We care about your success

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The sector of pharmacies, hospitals and medical practices is especially demanding in terms of temperature stability. Infrico Medcare in its Pharma CARE and Pharma CARE PLUS product lines has incorporated a series of specific characteristics that guarantee the safety in the conservation of medicines and vaccines 24 hours a day and 365 days a year. This is crucial for our customers.

All our refrigerators of the Pharma CARE PLUS and Pharma CARE PLUS lines are designed in compliance with DIN 58345.

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**According to DIN 58345 the devices must comply with the following requirements:**

- Operating temperature range from +2°C to +8°C.
- Interior room temperature of the installation site is accepted between +10°C and +35°C.
- Interior is easy to clean and disinfect; water resistant and with corrosion protection.
- Mechanical capacity load of uniformly mounted parts of 100 kg/m².
- Doors fitted with locks.
- Acoustic emission below 60 dB [A]
- Power failure alarm for a minimum of 12 h.
- Optical-acoustic alarm in case of temperature deviation.
- Potential free alarm contact.
- Safety device to prevent the temperature going below +2°C

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Using INVERTER technology in our equipment, we achieve greater temperature stability, lower noise level and outstanding energy efficiency.

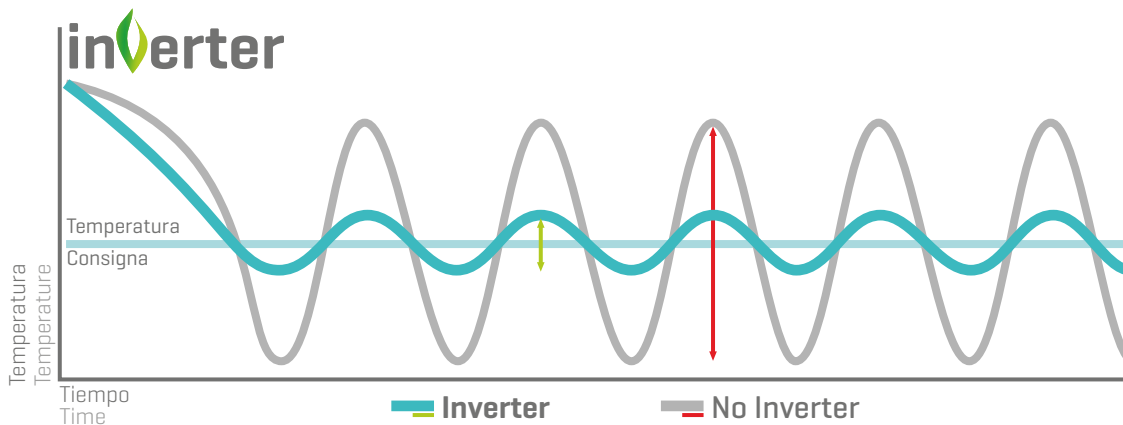
**MENOR CONSUMO ENERGÉTICO**  
**LOWER ENERGY CONSUMPTION**

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
Conventional non-INVERTER equipment, the compressor operates at maximum speed until reaching the setpoint temperature, and once it is reached, it stops, repeating successively with continuous starts and stops, causing greater energy consumption and lower temperature stability.

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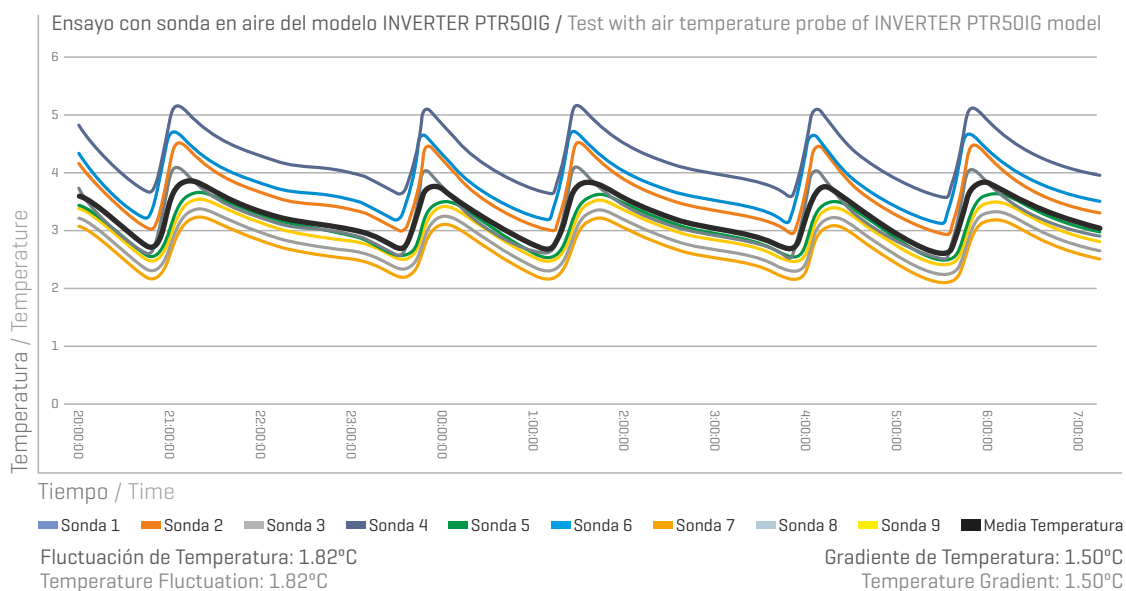In equipment with INVERTER technology, once the compressor starts operating, it adjusts the operational speed to the setpoint temperature, avoiding the starts and stops of the compressor that increase energy consumption and make greater the oscillation of the equipment temperature.

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OUR EXTENSIVE EXPERIENCE AS A WORLDWIDE MANUFACTURER OF COOLING EQUIPMENT AND THE SPECIFICATIONS OR QUALITY REQUIREMENTS OF DIFFERENT MARKETS THAT WE COMPLY WITH, ENSURES YOU CAN RELY ON INFRICO MEDCARE.

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Pharma CARE SERIES offers pharmaceutical refrigerators in different sizes and storage capacities, in order to fit all possible space requirements and needs. Models range from countertop, underbench or upright versions which are available with model variety from 70 to 360 litres.

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These models are designed according to **DIN 58435** standard.

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This series offers high performance refrigerators for the storage of pharmaceutical preparations to meet the most demanding applications. In order to achieve that purpose, these models have incorporated **INVERTER** technology with high energy efficiency and low energy consumption and lower noise level.

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
These models are designed according to **DIN 58435** standard. Capacity ranges from 170 to 1300 litres, and all models with Digital Plus. controller.

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Solid or glass doors are available in the whole range.

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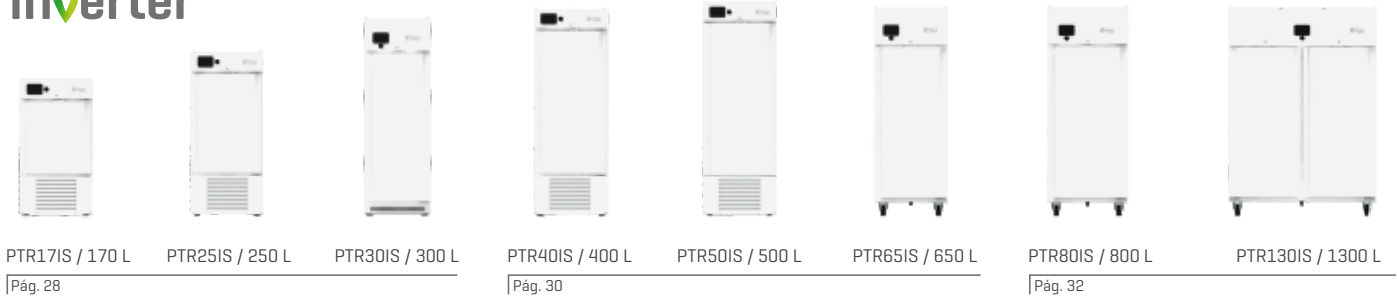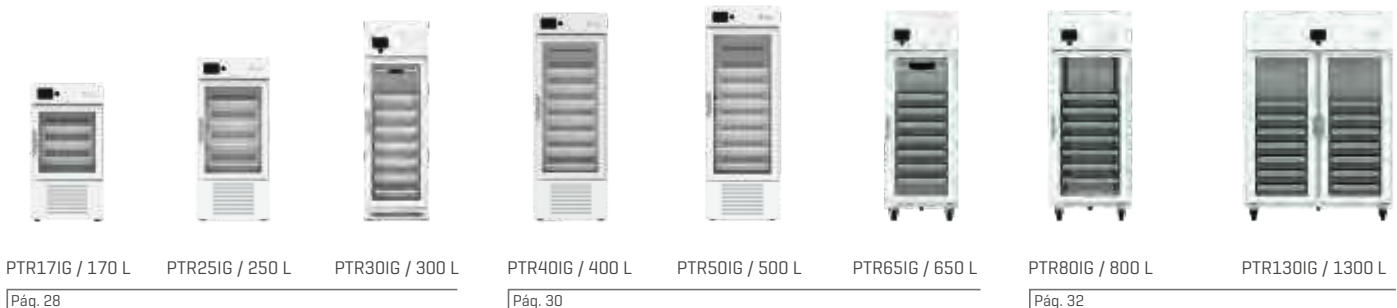
**DIGITAL CONTROLLER** (Pharma CARE range models] or **DIGITAL PLUS CONTROLLER** (Pharma CARE and Pharma CARE PLUS range models)  
Choose according to your needs!

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**LOCK / DOOR SENSOR**  
All our equipment has locks and open door sensors, to guarantee the security of the stored products by controlling opening time and fans functioning.

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**LED LIGHTING**  
The use of LED lighting reduces the energy consumption and improves the homogeneity of the temperature inside the cabinet.

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**FLEXIBLE STORAGE SOLUTIONS**  
Our new shelf / drawer system allows you to adjust their position, vary their number, change shelves for drawers and vice versa or compliment them with dividers. This provides the fit of the equipment to your needs and storage convenience of medications or vaccines of different sizes.

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**DOOR STYLE**  
It is possible to choose solid or glass doors in our whole range. All our doors are hinged and reversible.

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**EASY REMOVABLE DOOR GASKET**  
It is ideal to maintain optimal hygiene and facilitate maintenance work.

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**ROUNDED CORNERS**  
The interiors have rounded corners, which facilitate their cleaning.

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### BATTERY

Allows the maintenance of temperature and alarms register in case of a power failure.

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### INVERTER

INVERTER technology (Pharma CARE PLUS range models) with high energy efficiency and low energy consumption.

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### FREEZING PROTECTION ELECTRONIC CONTROLLER

Safety thermostat prevents the temperature dropping below +2°C.

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### EVAPORATION PAN

Our equipment has a built-in tray to evaporate automatically defrosted water.

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### ADVANCED DEFROST SYSTEM

Built-in to ensure uniform cooling over time and to facilitate maintenance. In addition, the equipment analyses its operation and only defrosts when it is necessary, to improve the stability of the temperature.

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### CASTERS / ADJUSTABLE LEGS / HEIGHT REGULATORS

In our whole range, there is a possibility to incorporate casters to facilitate displacement and cleaning.

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
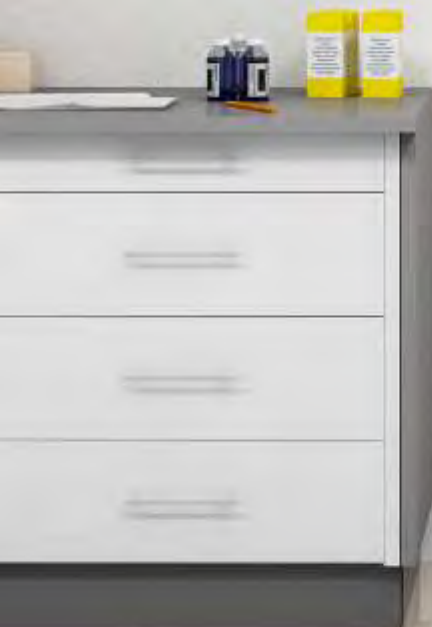
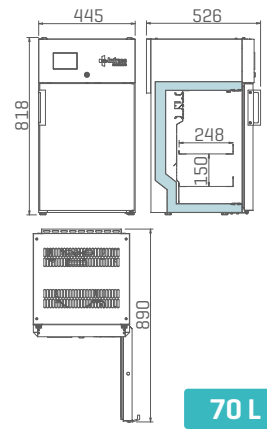
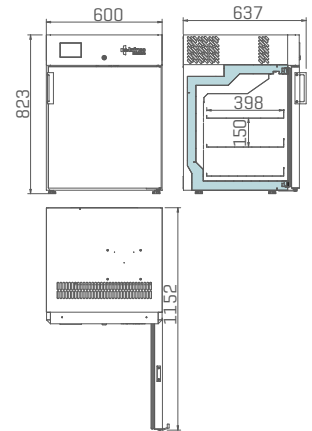
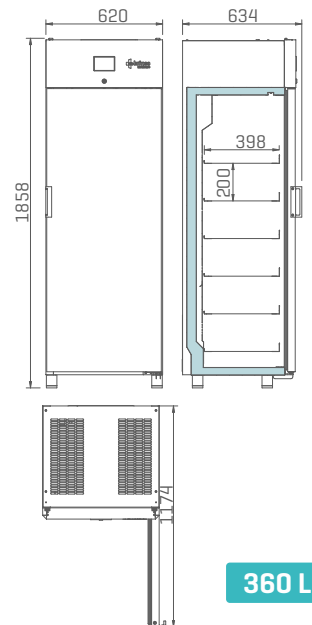
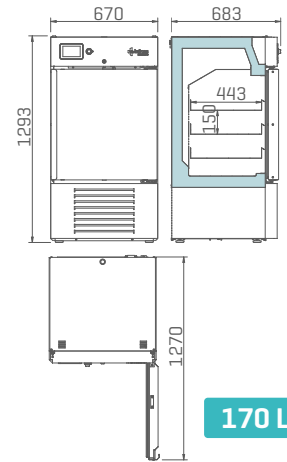
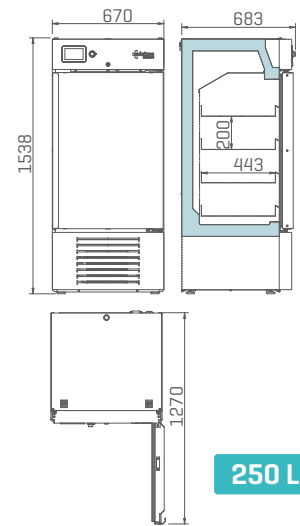
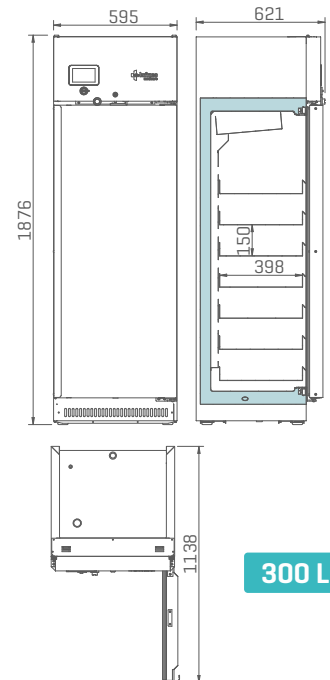
We reserve the right to modify technical and construction specifications without prior notice.

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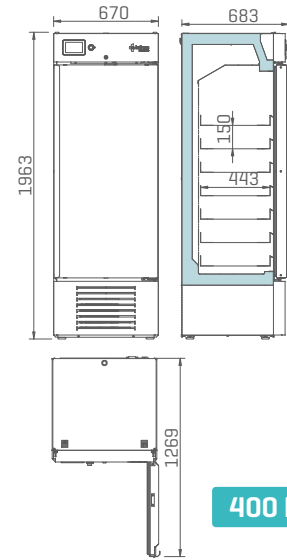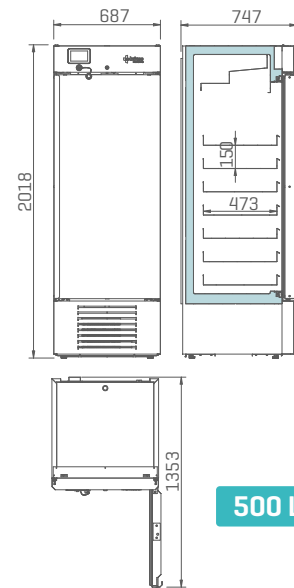
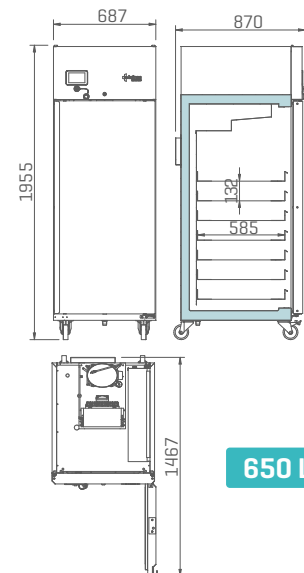
\* Measured at +25°C ambient temperature and

GAMA SERIES	Pharma CARE PLUS	Pharma CARE PLUS	Pharma CARE PLUS	Pharma CARE PLUS
MODELO MODEL	PTR80IS	PTR80IG	PTR130IS	PTR130IG
Rango T° Temperature range [°C]	+2 / +8	+2 / +8	+2 / +8	+2 / +8
Capacidad Nominal (litros) Nominal Capacity [litres]	800	800	1300	1300
Medidas Exterior [mm] L x P x A Exterior dimensions [mm] W x D x H	797 x 1000 x 1955	797 x 1000 x 1955	1385 x 834 x 1955	1385 x 834 x 1955
Medidas Interior [mm] L x P x A Interior dimensions [mm] W x D x H	645 x 803 x 1388	645 x 823 x 1388	1234 x 681 x 1388	1234 x 702 x 1388
Claro de puerta [mm] L x A Door opening [mm] W x H	645 x 1350	645 x 1350	534 x 1350	534 x 1350
Espesor Aislamiento [mm] Insulation thickness [mm]	75			
Acabado Exterior Exterior finishing	Acero galvanizado plastificado blanco White plastic coated galvanised steel			
Acabado Interior Interior finishing	Acero Inoxidable Stainless steel			
N° Cajones Drawers Number	7	7	14	14
Dimensiones Cajón [mm] L x P Drawer Size [mm] W x D	564 x 750	564 x 750	452 x 585	452 x 585
Material Cajones Drawers Material	Acero Inoxidable Stainless steel			
Intercambiabilidad de Estantes/Cajones Interchangeable Shelves/Drawers	Si Yes			
Tipo de Puerta Door Type	Ciega Solid	Cristal Glass	Ciega Solid	Cristal Glass
Cerradura Lock	Si Yes			
Salida USB USB Port	Si Yes			
Datalogger Datalogger	Si Yes			
Batería Battery	Si Yes			
Sonda simulación producto Product temperature simulation probe	Opcional Optional			
Sonda de temperatura ambiente Interior room temperature probe	Si Yes			
Termostato anticongelación Freezing protection controller	Si Yes			
Tipo de Controlador Controller type	Digital Plus			
Tipo de Refrigerante Refrigerant type	R290			
Potencia Eléctrica (W) Electric power (W)	434	442	471	510
Tensión/Frecuencia [V / Hz] Voltage/Frequency [V / Hz]	230 / 50			
Peso [kg] Weight [kg]	172	178	208	225

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